ibérica

Industrial laborum
lab integrated systems

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EN

PHYTO-PHARMACEUTICALS
PESTICIDES
OTHER PRODUCTS FOR
AGRICULTURE USE

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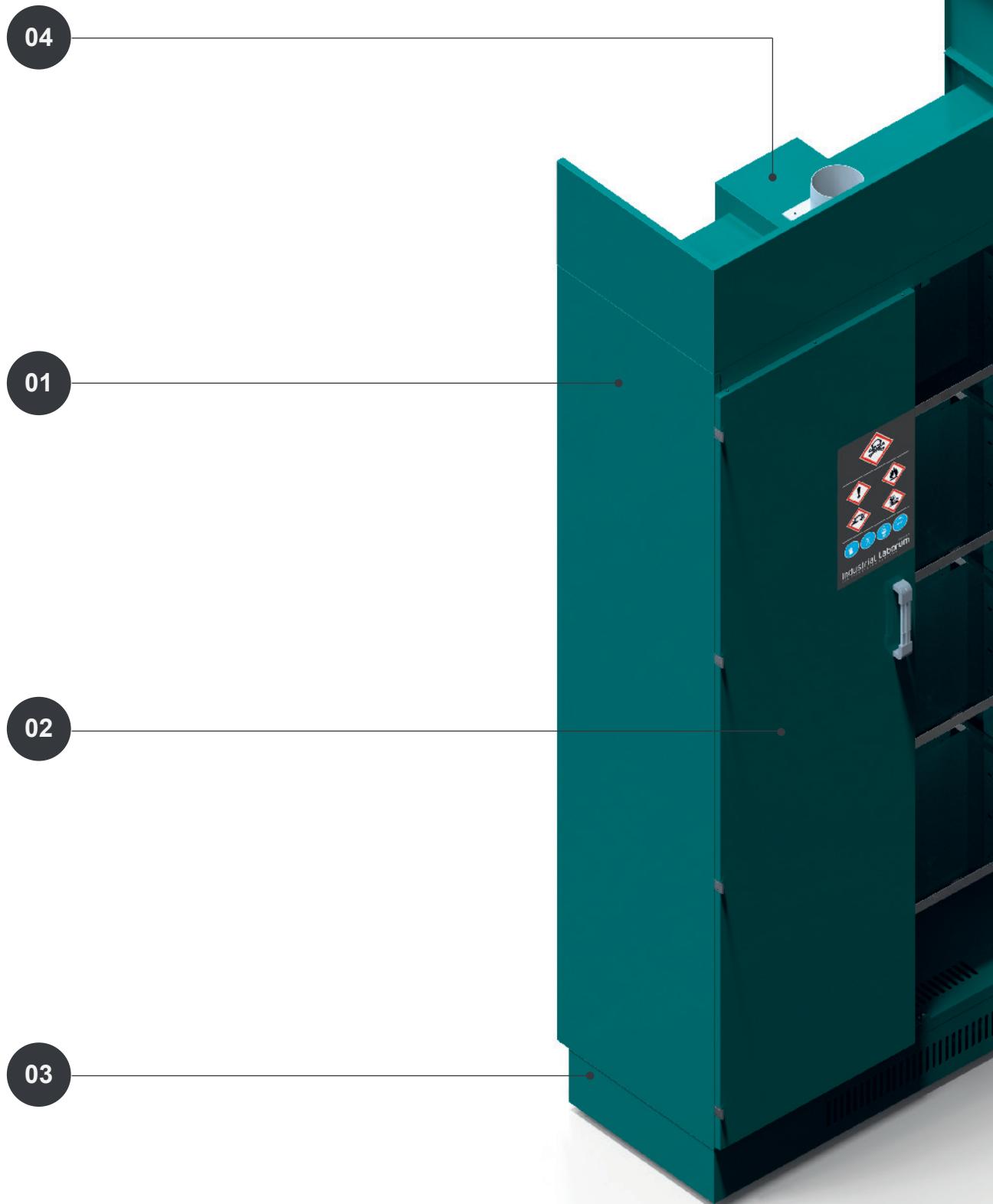
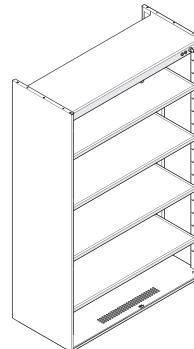
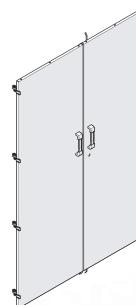
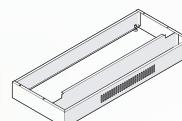
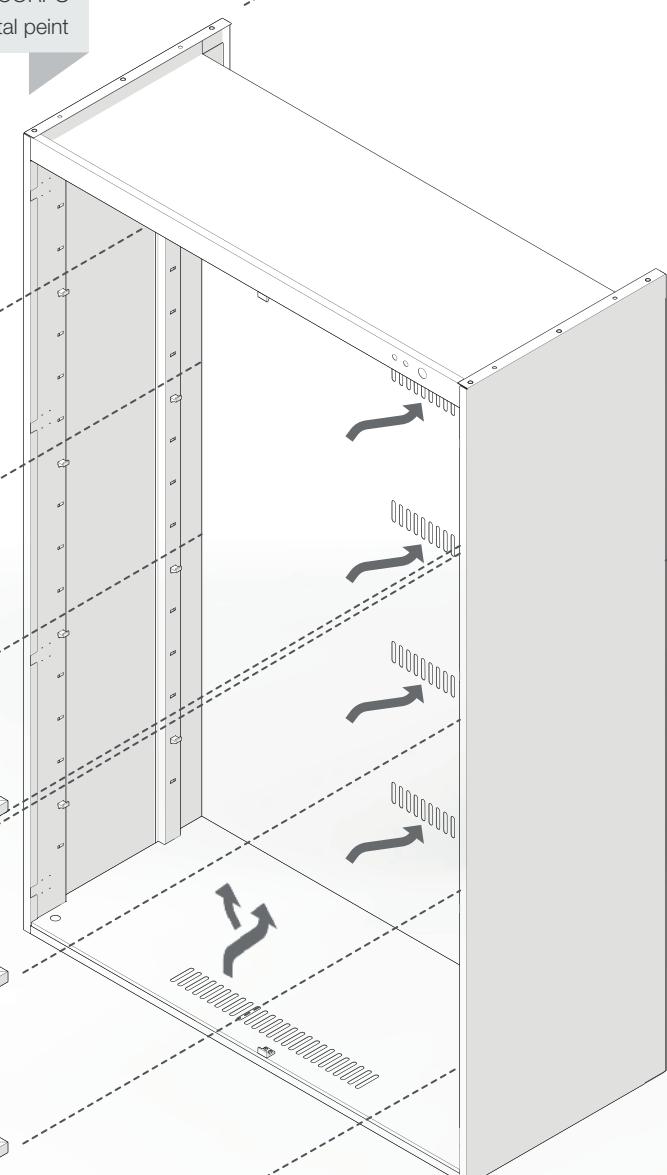
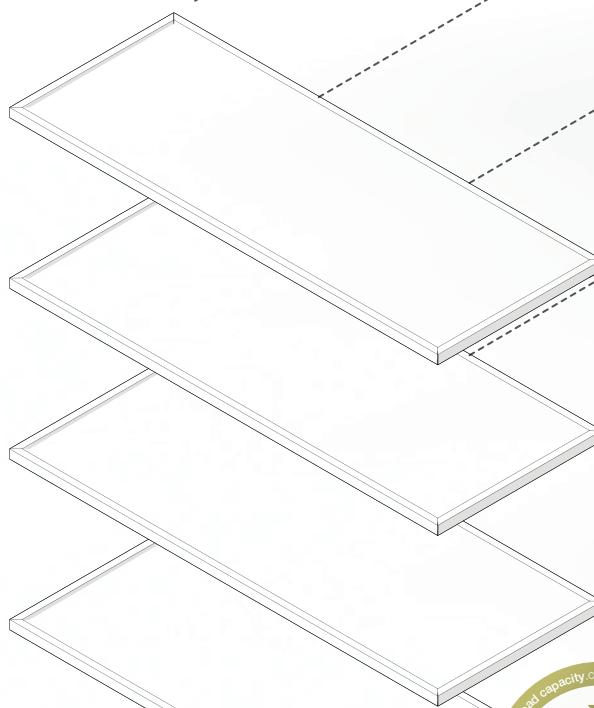
EN

CARCASE

The **carcasse** is welded constructed and manufactured with cold rolled steel sheet coated with thermosetting powder of epoxy-polyester resins.

The goal of the **openings in the bottom** is to allow the **entry of air** while the **back openings** are used to **exhaust the interior compartment** through the airtight air chamber, installed on the back of the cabinet.

The cabinet is equipped with 4 removable and height adjustable stainless steel **shelves**. All of them have **perimeter edge**, essential for the user's safety.

Maximum load capacity (evenly distributed): **15 kg** per shelf

Indoor and outdoor surfaces are smooth and uniform for an easy cleaning and disinfection.

High resistance to corrosion, abrasion and impact.

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160 mm

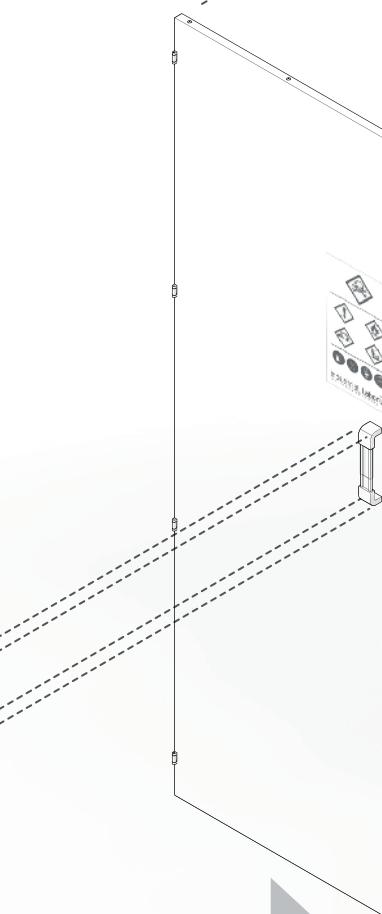
aluminium + plastique

DOOR
made out of painted metal
w/ sound absorbing filling

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EN

DOORS

Hinged opaque doors made out of cold rolled steel sheet panels, with expanded polystyrene **sound absorbing filling**, coated with thermosetting powder of epoxy-polyester resins.

Locking system with plastic components resistant to corrosion and bumpers for silent closure.

Smooth surfaces for easy cleaning and decontamination.

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The use of protective gloves is required.

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The use of protective clothing is required.

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The use of protective mask is required.

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TOXIC - May cause nausea, vomiting, headache, loss of consciousness or other damages, including death.

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IRRITANT [Xi] / HAZARDOUS [Xn] - May cause allergies, eczema, eyes, throat, nose or skin irritation. Exposure to high doses can cause somnolence or even poisoning.

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FLAMMABLE - May ignite in contact with flame, sparks, static electricity or heat exposure.

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CORROSIVE - Attacks or destroys the metals. May cause skin or eyes burns in case of contact or projection.

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INJURIOUS TO THE ENVIRONMENT - Toxic to aquatic organisms (fish, crustaceans or algae).

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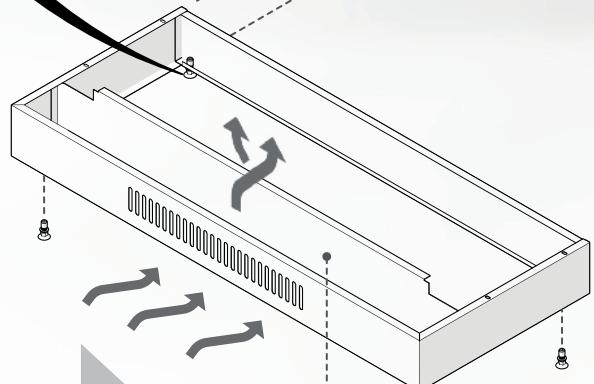
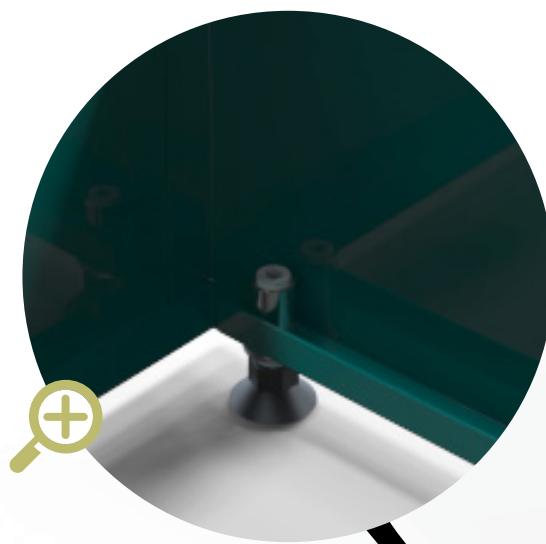
EN

PLINTH

Painted metal **closed plinth** with 4 polyamide **levellers** to compensate the floor irregularities.

The **openings** in the front allow **the entry of air** into the cabinet.

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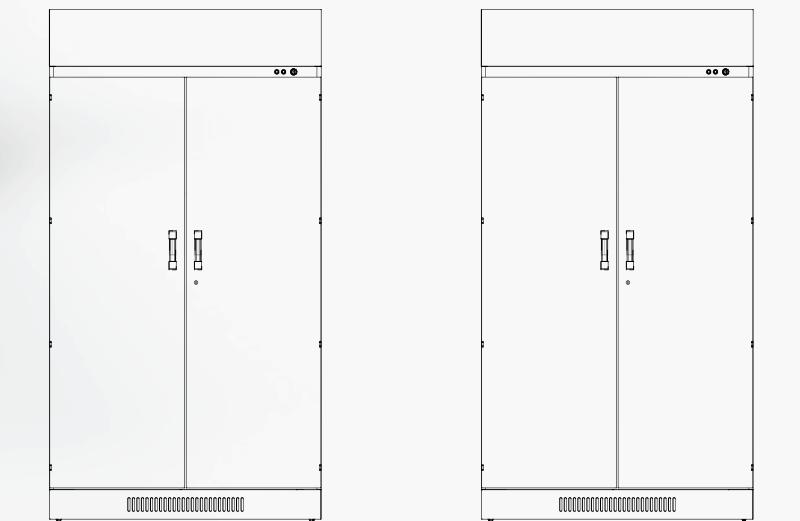
VENTILATION

The **cabinet** includes a **ventilation system** comprising:

- a painted metal sealed **collecting duct**, installed in the rear part of the cabinet, **to connect to the exhaustion**. It allows **the sweeping of the interior cabinet's compartments**;
- an **in-line ventilator** made out of polypropylene, **to be applied in the extraction outlet** of the cabinet;

The **ventilator's remote system** is composed by a **timer** (to automatically turn the ventilation on and off to ensure the minimum flow required and minimize energy consumption), a **control pushbutton (1)** and **two light signals on/off (2) (3)** installed in the upper part on the cabinet.

Once connected, the cabinet can **renew the inside air**, with the doors closed, **10 times per hour**.

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
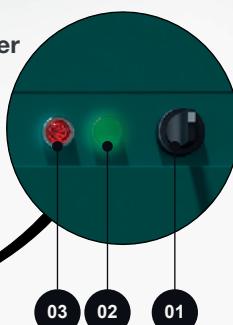
VENTILATION

EXTRACTION OUTLET
made out of polypropylene

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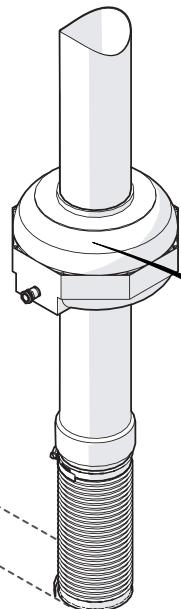FLEXIBLE SLEEVE + DUCTS
made out of polypropylene

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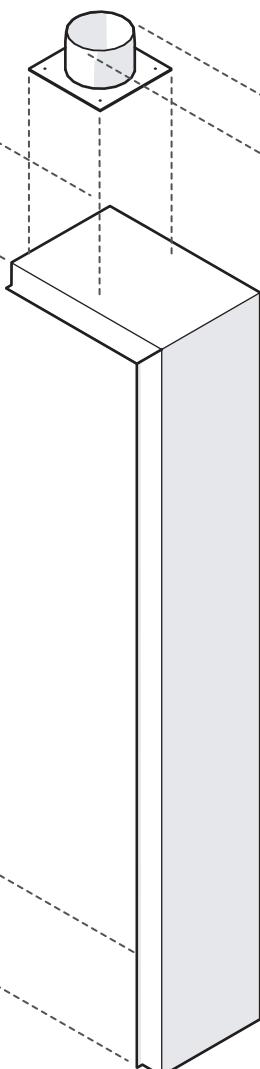
10 x 

CONNECTED CABINET
inside air renovation
10 times / hour

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
COLLECTING DUCT
made out of painted metal

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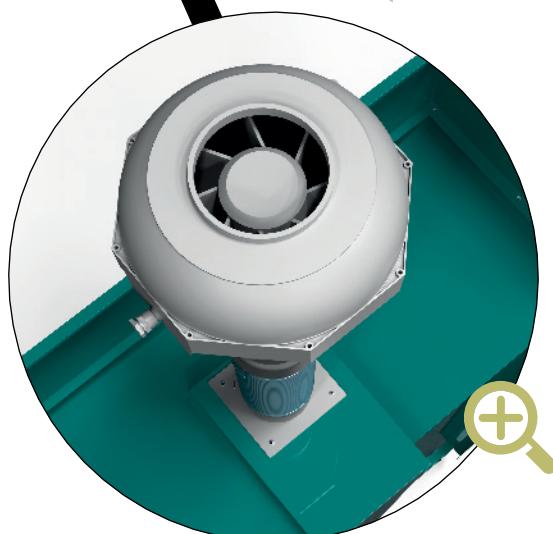IN-LINE VENTILATOR
made out of polypropylene

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EN

MATERIALS

PAINTED METAL EXTERNAL HOUSING *
STAINLESS STEEL SHELVES
PP VENTILATION SYSTEM

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